

# FLEXIBLE-ARM™ LIGHT INSTRUCTIONS

## SPECIFICATIONS:

Electrical: 150 VAC, 60Hz, UL Listed  
 Light: 35 watt Halogen bulb  
 Plug: Polarized  
 Arm length: **Cat. No. 3095** - 21"  
                   **Cat No. 3096** - 31"  
 Weight: 8 lbs.

## FEATURES

- Base - 7" diameter, 7lb.
- 35 watt bright Halogen white light
- Clamping unit for vertical or horizontal Positioning
- Switch in cord
- Flexible-Arm™ articulates at 3 points
- Fingertip bar for raising, lowering and pivoting with ease
- Spring counterbalance for stability
- Vented shade with rolled edges
- Safety-glass with UV filter

## BASE ASSEMBLY (see Figure 1)

1. Remove "Screw I" from base "Stem II".
2. Insert the lamp frame into base "Stem II", align screw holes.
3. Screw in "Screw I" and finger tighten.

## CLAMPING UNIT ASSEMBLY

1. Insert the lamp frame into hole in top of clamp assembly.
2. Loosen clamp with lever on bottom of assembly.
3. Position clamping mechanism and tighten firmly using lever on bottom of assembly.

## BULB REPLACEMENT (see Figure 2)

1. Turn off lamp and allow shade and accessories to cool.
2. Turn the lampshade upward.
3. Remove CENTER screw (using a Phillips screwdriver) and clip from protective glass cover.
4. Pull out failed bulb and replace with halogen bulb. (Control Company Cat. No. 3098)
5. Replace protective glass and clip and screw in place.

**NOTE:** Do not remove screws holding fingertip bar in place.

## WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

## CONTROL COMPANY

4455 Rex Road  
 Friendswood, Texas 77546 USA  
 Ph. 281-482-1714  
 Fax 281-482-9448  
 E-mail sales@control3.com  
 www.control3.com

Control Company is  
 ISO 9001 Quality-Certified by DNV and  
 ISO 17025 accredited as a  
 Calibration Laboratory by A2LA.

**Cat. No. 3095/3096**  
**Flexible-Arm™ Light Instructions**  
**92-3095-10**  
 Rev. 1 071504

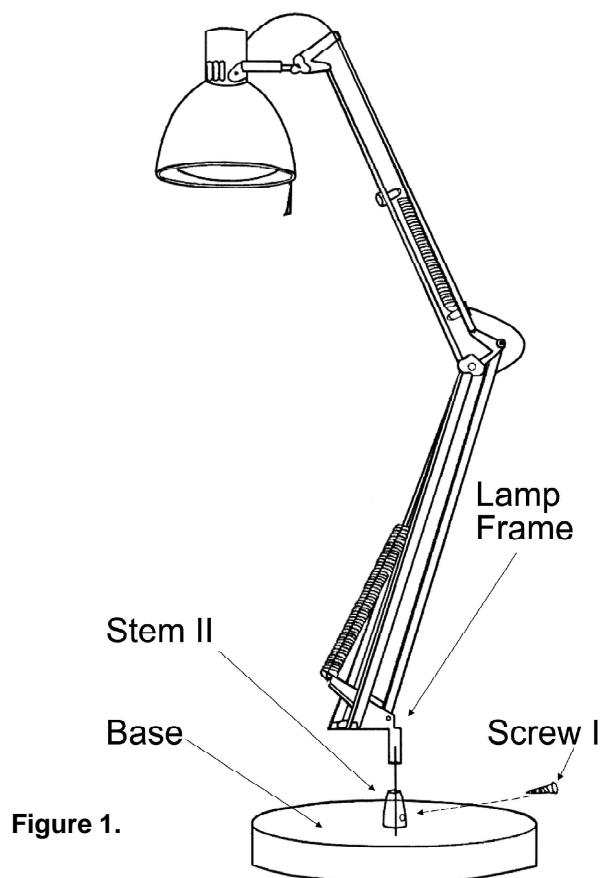


Figure 1.

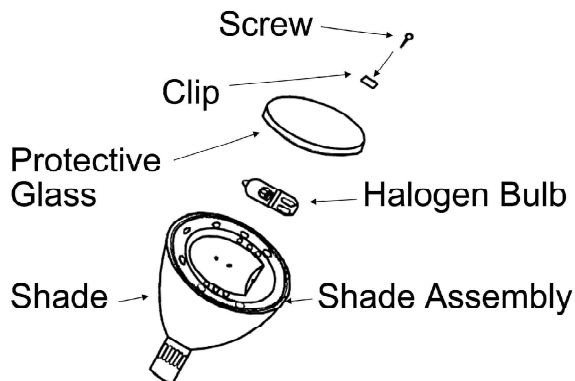


Figure 2.