

GLASSWARE RESCUER™ INSTRUCTIONS

SPECIFICATIONS

File length: 4 inches
File width: 1/4" half round
Handle length: 4 inches
Length of entire unit: 8 inches

Glassware Rescuer™ is constructed using real diamonds; the hardest material known. Use it to repair all beakers, flasks, and lab glassware, as well as most hard materials.

OPERATIONS

1. Moisten Glassware Rescuer™ with water.
2. With round side down, rub back and forth in the recessed edges of the broken object.
3. With flat side down, rub back and forth on the sharp raised corners of the broken object.

NOTE:

Only a few strokes should be needed to smooth most surfaces.

CAUTION:

Eye protection and gloves should be worn. Caution should be used when repairing all glassware.

CONTROL COMPANY

4455 Rex Road
Friendswood, Texas 77546 USA
Ph. 281-482-1714
Fax 281-482-9448
E-mail sales@control3.com
www.control3.com

Cat. No. 3315
Glassware Rescuer™ Instructions
92-3315-00
Rev. 1071707